

HAAKE Viscotester iQ OSC Update

The HAAKE Viscotester iQ rheometer is equipped with an optional oscillation mode which can be activated any time by applying the OSC Update (order number 222-2207).

Note The HAAKE Viscotester iQ must be equipped with firmware versions V1:01.07.000 and V2:1.04.000 or newer.


Applying the OSC Update

The OSC Update consists of an installation key which must be entered in a dialog using the HAAKE Viscotester iQ Update Tool.


See the “HAAKE VTiQ Reference Manual, Appendix D, Firmware update”, for detailed instructions on where to find, how to start and how to use the HAAKE Viscotester iQ Update Tool.

❖ To enter the installation key


1. Launch the HAAKE Viscotester iQ Update Tool (version 1.3.0.0) or newer.)

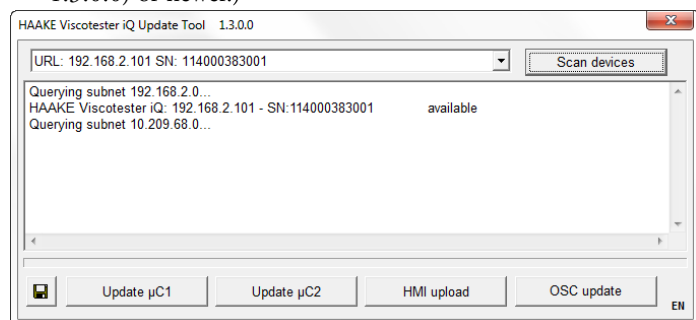

2. Click the **OSC update** button to open the OSC Update dialog. The number behind Connected is the serial number of the Viscotester iQ which must match the serial number in the installation key. The text OSC OFF indicates that the oscillation option is not yet activated.

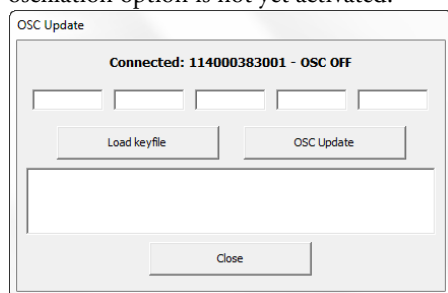


3. From the information on the sticker below, enter the key information in the five edit fields above the two buttons.

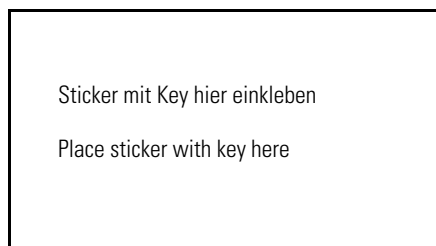

or


Click the **Load keyfile** button to read the key information from a key file if available.

4. When the update was successful the text SUCCESS OSC update will be displayed in the area below the two buttons.

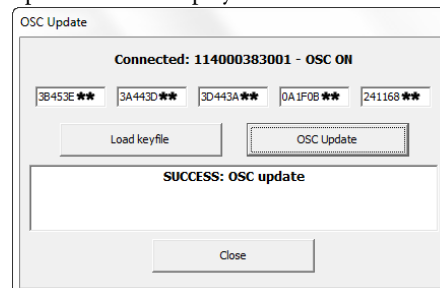

5. Click the **Close** button to close the OSC Update dialog.

A confirmation text for the update is now displayed in the HAAKE Viscotester iQ Update Tool program window.

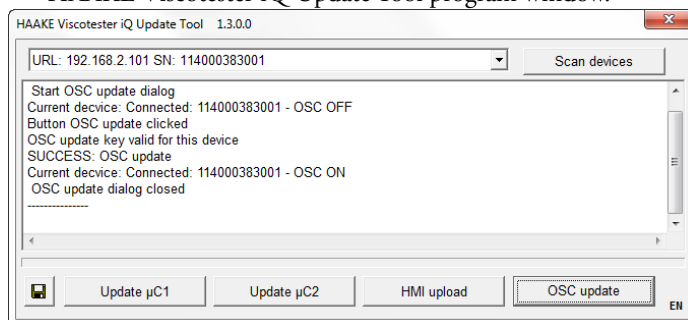

6. Click the **Close** button , to close the HAAKE Viscotester iQ Update Tool program window.

7. On the Viscotester iQ touchscreen panel tap the **Configuration** button.

8. In the Configuration menu select **Device info**.

In the first line of the Device info menu the text OSC should now be displayed to show that the oscillation option is activated.



Contents of delivery

Part	Number
HAAKE Viscotester iQ OSC update installation key	098-5073
This document	006-2764