

What's new in RheoWin

Version 4.95.0003

(Released 22.07.2025)

Below is a list of things that are new and/or changed in version 4.95.0002 compared with version 4.95.0000 of RheoWin.

Installation

Installation of iQUpdate.exe

- In RheoWin 4.95.0000 the iQUpdate.exe program which is needed to update the VT iQ (Air) and MARS iQ (Air) firmware was missing. This has been solved.

Measuring geometries

- Certain measuring pressure cell measuring geometries and the adapters PC1 and PC2 were not installed correctly. This has been solved.

DeviceManager / Device Drivers

Measuring Geometries

- The following pressure cell geometries are now part of the installation CC20-PC/SS, CC25-PC/Ti/SE, CC31,9-PC/Ti and CC38 DG-PC/Ha.
- All already available pressure cell geometries are now also listed with their new names which use the format CCxx-PC/Ti, CCxx-PC/Ha, CCxx DG-PC/Ha, FLxx yy-PC/Ti etc, where xx stands for the rotor diameter and yy for the vane blade height. The letters PC in the new names stands for Pressure Cell.

JobManager

All OSC Elements

- In RheoWin 4.95.0000 in all OSC elements the setting **Auto: High quality acq. (f-dependent)** for the **Waiting time**, **#Repetitions** on the **Data acquisition** page of the element editor, did not work correctly with the MARS 40/60 (and older MARS I/II/III). This has been solved.